T° SENTRY MODEL 140 Digital Alarm Module

OPERATING INSTRUCTIONS MANUAL

TIP TEMPerature Products

340 W Broad Street Burlington, NJ 08016 1-800-TIP-TEMP www.tiptemp.com

TABLE OF	CONTENTS
----------	----------

Introduction	1
UNIT CHECK	2
PANEL DESCRIPTION	3
UNPACKING INFORMATION	5
OPERATIONAL CHECK	5
BASIC OPERATION	6
SENSOR PROBE PROBE INSTALLATION PROBE LOCATION PROBE CALIBRATION	7 7 8 9
OPTIONAL FEATURES FAHRENHEIT CALIBRATION AND DISPLAY (F) HIGH RESOLUTION DISPLAY (H) RELAY (R) NIMH BATTERY BACK-UP (B) CURRENT OUTPUT FOR RECORDER (I) VOLTAGE OUTPUT FOR RECORDER (V) ANALOG COMPUTER MONITORING SYSTEM (CMS) DIGITAL COMPUTER MONITORING SYSTEM (D)	12 12 12 13 13 14 14
PROGRAMMING LIST OF PARAMETERS	15 16

LIMITED WARRANTY

TIP TEMPerature Products warrants each manufactured item against defects in material and workmanship, when used as recommended, for a period of one year from original purchase. Products believed to have such defects must be returned to the factory by prepaid transportation.

TIP TEMPerature's obligation under this warranty is limited to the repair or replacement, at its option, of those items which upon examination prove to be defective. Such repair or replacement will be made without charge.

This warranty will be void if repairs or alterations are made or attempted without factory authorization; or if the item has been subject to misuse, negligence or accident.

TIP TEMPerature Products assumes no liability for consequential damages of any kind. The purchaser, by acceptance of the product, assumes all liability of the consequence of its use or misuse.

TIP TEMPerature Products makes no other warranty, whether expressed or implied, in connection with the sale or use of its products.

INTRODUCTION

The Model 140 Digital Alarm is a temperature monitoring and alarm device with user-programmable high/low alarm set points and initial alarm delay time. Parameters are stored in non-volatile memory, and are maintained even when power is removed.

The temperature readout for the Model 140 can factory-set to °C or °F. Its small size makes it easily interchangeable with the Hampshire Controls Model 125 Digital Alarm Module. It can also be readily mounted into optional custom panels, for integration with existing equipment.

The Model 140 is very simple to operate: plug in the unit, place the probe in the location to be monitored, and the display will show the temperature of that location. Pushing three buttons allows the user to program both high and low alarm set points for the temperature range.

Optional features allow the unit to send temperature data to a recorder or (through an appropriate translator), a computerized monitoring system. Please contact Hampshire Controls for details of the many possibilities!

UNIT CHECK

Check the serial number label on the rear of the unit to be certain that it is the proper unit for your application. The label will show model number, temperature range, probe type, and any options in the following format: 140-T-P-O.

Temperature ranges:


T Code	Range, °C	Range °F
100	-100 to +100 °C	-148 to 212 °F
200	-200 to +50 °C	-328 to 106 °F
400	0 to +400 °C	32 to 752 °F

Probe Types:

PC	ode	Probe
P1	00	100Ω RTD (standard probe)
P1	000	1000Ω RTD
- 1	J	Micro-amp current source
1	K	K-type thermocouple
	Т	T-type thermocouple

Options:

Option
Fahrenheit display option
High resolution display (0.1°) (not all ranges or sensors)
Relay output
NiMH battery back-up with built-in trickle charger
4-20 ma current output
Voltage output for recorder
Analog computer monitoring system (used w/CMS)
Digital serial communication
RS232 or RS485 or CMS translator
Custom option

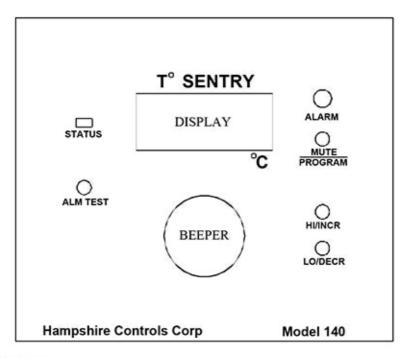

PANEL

The <u>LCD Display</u> shows the temperature in °C or °F, depending on the option.

 If the unit has the battery-backup option, the display shows a BAT message to indicate that there has been a power loss and it is running on the battery.

The STATUS tricolor LED shows the status of the unit.

Green = temperature in range

Red = temperature is out of range

Amber = program mode

The light blinks off briefly each second to show the unit is active.

The <u>ALM TEST</u> push-button interrupts line power, to allow testing of the battery and the alarm relay.

The bright red ALARM LED shows alarm conditions.

- In alarm mode, it flashes every half-second.
- In alarm mute mode, it gives a short blink every second.

The BEEPER shows alarm conditions and user activity.

- In alarm mode, it gives a long beep.
- In alarm mute mode, it gives a short chirp.
- It gives an audible tick to indicate a button has been pressed.

The MUTE/PROGRAM button has three functions.

- In Normal Run mode, if the unit is beeping, pressing the button will mute the beeper for a preset time. Each time the button is pressed the mute time resets.
- Holding the button down for two seconds takes the unit out of Normal Run mode and into Programming mode.
 At the same time, it mutes the beeper.
- When in Programming mode, pressing the button steps the unit through the various programming parameters.

The <u>HI/INCR</u> button has a different function for each operating mode.

- In Normal Run mode, pressing the button causes the high temperature set point to be displayed.
- In Programming mode, pressing the button increases the displayed parameter value.

The <u>LO/DECR</u> button has a different function for each operating mode.

- In Normal Run mode, pressing the button causes the low temperature set point to be displayed.
- In Programming mode, pressing the button decreases the displayed parameter value.

T° SENTRY MODEL 140 OPERATING INSTRUCTIONS

UNPACKING INFORMATION

The 140 unit normally ships with the probe connected.

If the unit has the battery back-up option, the battery is disabled before shipping.

NOTE: Remove the tagged screw in the side of the 140 box to enable the battery. Assume the battery is discharged, and allow it to charge for at least 24 hours to ensure full capacity.

OPERATIONAL CHECK

- Plug the Model 140 into a 110-120 volt power outlet. If the air temperature is within the unit's operating range, the display should show the current probe (air) temperature.
 - If the display is blank, check to make sure that power is going to the unit. If the STATUS LED is lit (any color), the Model 140 has power. If not, verify that the power supply is plugged into the unit and into line power.
- Hold the sensing part of the probe tightly in your hand. If your body temperature is within the unit's operating range, the displayed temperature should begin to go up. It will take several seconds to level off.
- If the Model 140 does not respond as it should, call Hampshire Controls for assistance.

BASIC OPERATION

The Model 140 has two modes: operating and programming. This section describes the unit in Operating Mode.

In operating mode, the display shows the temperature that the probe is reading. The display does not "round off" the temperature. If it shows a value of 10°, the actual temperature can range from 10.0° to 10.9°. Therefore, a very small system change of 0.1° or less (or a minor electronic disturbance) may cause the displayed temperature to change one digit. It is only when the displayed change holds steady for more than 20 seconds that it can be related to real thermal changes in the probe.

Press the push button labeled "HI/INCR" to cause the display to show the high temperature set point.

Press the push-button labeled "LO/DECR" to cause the display to show the low temperature set point.

Press the MUTE and DECR buttons at the same time to cause the display to show the software version number. It will be different from the temperature display in that it will show two decimal places, such as "1.01".

If the unit has the "Switch" option, press the HI/INCR and LO/DECR buttons together to cause the display to show if the switch is open ("OP") or closed ("CL"). When the buttons are released, the display will go back to showing the temperature.

The Model 140 compares the probe's temperature reading to the set points for out-of-range states. An out-of-range state is a temperature that is "equal to or beyond the set point". For instance, if the low set point is 0.0, then the temperature reading is out-of-range for 0.0° and below. The STATUS LED glows green when the probe reads an in-range temperature and red when the temperature is out of range.

T° SENTRY MODEL 140 OPERATING INSTRUCTIONS

The Model 140 will not declare an alarm until the out-ofrange state has continued for a specified time (the "alarm delay"). This alarm delay time will prevent transient nuisance alarms.

When an alarm condition occurs, the beeper and LED will both begin. Press the MUTE/PROGRAM button to stop the beeper for a specified time, changing the sound to short 'chirps'. When the temperature goes back within range and there are no out-of-range readings for thirty seconds, the chirping stops and the MUTE/PROGRAM button resets. If another out-of-range condition then occurs, the beeper will again start at full volume.

After the user-programmable "SILENCE" parameter times out, if the Model 140 is still in a constant alarm condition, the beeper will go back to full volume.

SENSOR PROBE

The probe supplied with the Model 140 is a highly accurate temperature sensor. The sensor may be put in water or other mild aqueous solutions.

NOTE: Avoid submerging the probe in solvents or harsh chemicals. Use protective thermowells when monitoring such materials.

NOTE: The warranty does NOT cover damage to probes or electronics that is caused by exceeding temperature limitations, or damage to probes caused by using them in solvents or other unsuitable environments.

NOTE: If the display shows a continuous reading of 199.9, the probe has failed. Replace the probe and recalibrate the unit.

Probe Installation

The probe may be used in air or in simulated product.

If installing the sensor in a refrigerator, try to place the probe in a water and glycol chlorine solution to stop bacteria formation.

If installing the sensor in a cabinet or enclosure (particularly freezers), make sure to use good techniques to prevent room moisture from getting into the cabinet. Whenever possible, install the probe through an existing access port provided by the cabinet manufacturer, then reseal the port.

Alternatively, the probe can be run under, over, or through the door-sealing gasket. Often a door gasket will have a joint at one or more corners. Open that joint slightly by carefully making a slit with a razor blade. Insert the probe wire and then reseal the joint with flexible silicone sealing compound. Inside the cabinet, run the probe wire so that it will not become snagged during loading, unloading or cleaning procedures.

Probe Location

Install the sensor probe in a location where it will respond to the average temperature of the space being monitored and not to local conditions caused by door openings, etc.

The object of the probe location is to provide some safety for the area being monitored without generating "false" or nuisance alarms. For example, locating the sensor probe on the bottom of a chest freezer will result in the alarm being sounded later than if it was located near the top. However, locating the sensor too close to the top of the chest freezer could result in the alarm being sounded due to routine lid opening. Choose a probe location that offers the safety desired for the enclosure contents

Probe Calibration: Software Offset

Place the T140 probe and a reference probe together, in a stable well agitated bath. It is imperative that the reference probe and the T140 probe be positioned together and given time to equilibrate. If you find that the T140 temperature differs from your reference temperature, you may adjust a software parameter to offset the T140 temperature.

To access the OFFSET parameter

- Press and hold LO/DECR, then add HI/INCR, then add MUTE/PROGRAM,
- release LO/DECR
- Wait for the unit to give a short BLIP and the display to change to "OFF".
- Release the buttons.

Set the parameter accordingly. If you need the unit to display 2 degrees lower than it does, set the OFFSET to -2. If you need it to display 1 degree higher, set it to 1.

Press MUTE/PROGRAM to store the parameter and advance to the next parameter. Press MUTE/PROGRAM to step out of Edit mode and back to the Run mode.

Probe Calibration: Full Hardware Calibration

The temperature sensor used in the Model 140 has excellent long-term stability and should not need recalibration in normal usage. However, if the probe is subjected to temperature extremes outside of the normal operating range for the unit, or if the probe is damaged and needs to be replaced, it must be recalibrated.

To recalibrate the Model 140, follow these directions:

Equipment Needed:

- Small flat screwdriver or potentiometer adjuster
- Water bath, in an insulated container and at a KNOWN temperature different from 0 °C. Use a thermometer you can trust.

- Ice water bath, with plenty of clean solid ice in it. The following bath will provide a temperature of 0°C to within ±.04°:
- Use an insulated Thermos container to contain the bath.
- Get an ordinary aquarium bubbler. It should have a "stone" in it to make little bubbles rather than big ones. Put the end of the bubbler hose in the bottom of the Thermos container.
- Place cold water in the container, with plenty of ice. The water and ice should be distilled water, since water with impurities in it freezes at a temperature below 0°C.
- Start running the bubbler.
- Use this water bath rather than just putting the sensor on plain ice because the plain ice will probably be <u>colder</u> than 0°C (freezers generally operate at –20°C). The air from the bubbler keeps the water moving over the ice. The ice chills the water, but since the water is kept moving by the bubbler it doesn't freeze. The cold of the ice keeps it right at the freezing point, which for distilled water is 0°C.

1. Set up the unit for calibration:

- 1.1. Disconnect the unit from its power source.
- 1.2. Unscrew the four corner screws on the front panel. Carefully remove front panel from box. Turn it over and lay it down to the right of the case. The circuit board of unit is exposed to view.
- 1.3. Plug the unit into a 110-120 volt power supply.

2. Adjust the Zero Point:

2.1. Stir the ice water bath and place the temperature probe in it. Let the probe sit in the bath for about 5 minutes to allow it to stabilize at temperature.

- 2.1. Locate the adjustment potentiometers (pot) on the left edge of the board. One of the adjustment potentiometers will be marked with a "0". This is the ZERO adjust. Another will be marked with an "S". This is the SPAN adjust.
- 2.2. Turn the "0" adjustment pot with a small screwdriver until the display reads 0 (or 0.0 for H option models). Try to "center" 0 by turning a little past until a 1 or -1 appears, estimating the distance moved, then returning the pot to a point where 0 just appears on the display.

3. Adjust the Span:

- 3.1. Using a rubber band or masking tape, join the sensor probe and the thermometer so that their sensing tips are as close together as possible.
- Stir the probe/thermometer in the water bath. Allow time for the probe and the reference thermometer to stabilize.
- 3.3. Turn the "S" adjustment pot until the reference thermometer and the display agree. Use the same "centering" technique outlined under the zero adjust instructions.

NOTE: If this higher temperature is changing rapidly, it will be difficult to make an accurate adjustment. One technique to slow the rate of change is to use an insulated container (Dewar flask or Thermos bottle), and keep the probe/thermometer near the bottom while stirring gently.

4. Repeat 2 & 3

The SPAN and ZERO adjustments are interactive, therefore recheck the ZERO setting. Readjust if needed and then recheck the SPAN adjustment.

NOTE: Factory recalibration of units that have been incorrectly adjusted is not covered by warranty

OPTIONAL FEATURES

Fahrenheit Option (F)

The standard 140 displays temperature in degrees Celsius. The "F" option indicates that the unit is calibrated and displays degrees Fahrenheit.

High Resolution Display (H)

The high-resolution option allows temperature monitoring and alarm setting to 0.1°C. All of the Model 140 standard features remain the same.

Relay Option (R)

The relay option consists of a single-pole double-throw (SPDT) relay that changes state (switches) with the loss of line power or with the presence of an alarm condition.

NOTE: The relay is rated for pilot duty operation only. Do NOT use it for control switching. The relay contacts and connections are rated at 30v/1A.

In normal applications, the relay signals remote or central monitoring systems that an alarm condition has occurred at the equipment being monitored.

If the 140 unit has the relay option, the relay will "pull-in" when line power is present. If an alarm condition lasts for a designated time (the "relay delay"), the relay will "drop out", causing the relay contacts to switch. If the alarm condition does not last long enough for the relay delay to time out, the relay will not go into its alarm position. The relay will also "drop out" if power to the alarm is lost, acting as a power loss detector. When the main power to the alarm is lost, the bicolor LED on the upper left of the panel will go dark.

NOTE: Mute button, beeper or LED activities do not affect relay status.

Battery Back-Up (B)

Model 140 units that have the battery back-up option ship with the battery in place and connected, but disabled. Remove the screw located on the outside of the left side of the module to activate the battery option.

When line power is present, all of the normal features of the Model 140 are operating. When line power is lost, the battery provides power to continue to display the temperature of the area being monitored.

Testing the battery:

The battery recharges when the unit is plugged into line power. To test the battery, push the ALM TEST button on the front panel. The following should happen:

- The display will show a small BAT designation.
- The Alarm LED will light red.
- The beeper alarm will sound full on.
- On units with the relay option, the relay will switch.
 If the button is held longer than 6 seconds, the unit will reset,
 the activity LED will stop blinking, and the unit will not restart
 until 5 seconds after the button is released.

Current Output Option (I)

The I option sends out a milliamp signal that is proportional to the temperature being monitored. The I option is calibrated so that the units sends out 4 ma at a user defined low temperature, and 20 ma at a user defined upper temperature.

Voltage Output Option (V)

The V option sends out a millivolt signal that is proportional to the temperature being monitored. The standard output is 10 mv per degree C, with 0 mv at 0 °C. For example, if the monitored temperature is +37 °C, the voltage output is +370 mv (0.370 v). If the monitored temperature is -20 °C, the voltage present at the V option connector is -200 mv (-0.200v).

NOTE: This voltage output signal is intended for recorders and data loggers only. Do not try to use this voltage to operate other devices directly.

The input impedance of the apparatus using the voltage signal should not be less than 100K ohms.

Analog Computer Monitoring System Option (CMS)

This computer monitoring option is used with our computer based CMS system. This option provides a 4 prong terminal strip for power in and voltage out.

<u>Digital Computer Monitoring System (D)</u>

This serial communication option can be configured for RS232, RS485, or RS485/DMS. Please contact Hampshire Controls for serial protocol.

PROGRAMMING

To program the Model 140:

- Hold the Mute button for two seconds to bring the unit into Programming mode.
- The display will show the first parameter that can be programmed (high temperature limit). The name of the parameter will flash, then the current value of that parameter.
- To change the value of a parameter, press the INCR or DECR buttons. Pressing either button will change the parameter value by one count.
- Holding either button for half a second will change the parameter value by 10 counts. It will continue to step by ten on further button pushes, as long as the pushes occur within a short period.
- Releasing the INCR or DECR buttons for a half second will cause the step to revert back to one.
- 6. Once the unit is in Programming mode, each time the Mute button is pressed, the unit steps to the next parameter. As the unit steps to the next parameter, the value of the previous parameter is saved. Once the final parameter is reached, pressing the Mute button returns the unit to Normal Run mode, saving the final parameter.

Description of Parameters:

Standard Parameters:

Hi, Lo, Ad, rd, SIL

Hi (High Limit)

This is the high temperature limit. The current temperature is considered out-of-range if the probe reads a temperature equal to or above the high temperature limit.

Lo (Low Limit)

This is the low temperature limit. The current temperature is considered out-of-range if the probe reads a temperature equal to or below the low temperature limit.

Ad (Alarm Delay)

The Model 140 will not declare an alarm until the out-ofrange state has continued for a specified time (the "alarm delay"). This alarm delay time will prevent transient nuisance alarms. The alarm delay time can be set for 0 to 30 minutes.

rd (Relay Delay)

This parameter is used in conjunction with the relay option. If there is an alarm condition, the relay contacts will switch. The relay delay parameter allows the user to set the amount of time delay from when the alarm condition begins to when the relay contact actually switches. This keeps the relay from switching during short "nuisance alarms". The operator may set the relay delay time for 0 to 30 minutes.

SIL (Alarm Silence (Mute))

This parameter sets the number of minutes the beeper will "chirp" when the mute button is pressed during an alarm condition. After the Silence Time elapses, if the alarm condition still exists, the beeper will begin to sound at full volume again. This parameter can be set from 5 to 120 minutes.

Quick List of Button Functions

While in Run mode:

Press:	To:
HI/INCR	Display high temperature alarm point
LO/DECR	Display low temperature alarm point
MUTE & LO/DECR	Display software version number
HI/INCR & LO/DECR	For units with "Switch" option, display whether switch is open ("OP") or closed ("CL").
MUTE/PROGRAM	Mute alarm beeper for a short time.
MUTE/PROGRAM for 2 seconds	Bring unit into Programming mode for standard parameters
MUTE/PROGRAM & HI/INCR for 2 seconds	Bring unit into Programming mode for setup/special parameters only.

While in Programming mode

Press:	To:
HI/INCR	Increase parameter value
LO/DECR	Decrease parameter value
MUTE	Advance to next parameter